

# High Tryptone Cytophaga (HTC)

## Détails

For 1 liter of HTC:

4 g of Tryptone.

0.5 g of Yeast extract.

0.5 g of Beef extract.

0.2 g of Sodium acetate.

Complete to 1 liter with water.

pH7.3.

Autoclave.

Add 3 mM of sterile  $\text{MgCl}_2 \cdot 6\text{H}_2\text{O}$  and 10 mM of sterile  $\text{CaCl}_2$ .

For HTC solid medium:

Add 5 g of gelrite per liter.

Last modifications: 2011-02-21