

Alcalophilus medium (pH10)

Détails

For 1 liter of broth:

Part A: 900 ml of water 10 g of KH_2PO_4 5 g of peptone 5 g of yeast extract 1 g of glucose 0.2 g of $\text{MgSO}_4 \cdot 7\text{H}_2\text{O}$	Part B: 100 ml of water 20 g of Na_2CO_3
---	---

For agar:

Add 15 g of agar per liter to the part A

For top agar:

Add 7.5 g of agar per liter to the part A

Autoclave.

Mix the part A and part B.

Slepecky, R. A., and H. E. Hemphill. 1992. The Genus *Bacillus* - nonmedical, p. 1663-1696. In A. B. e. al. (ed.), *The Prokaryotes*, 2nd ed. Springer-Verlag, New York Inc., Secaucus, New Jersey. Dernières modifications: 16 janvier 2011