

## Media recipe

# TYA

### Détails

For 1 liter of TYA:

6 g of Tryptone.

2 g of Yeast extract.

3 g of Ammonium acetate.

0.2 g of  $\text{MgSO}_4 \cdot 7\text{H}_2\text{O}$ .

0.01 g of  $\text{FeSO}_4 \cdot 7\text{H}_2\text{O}$ .

Complete to 1 liter with water.

Autoclave.

Add 1% of sterile glucose and 3.7 mM of  $\text{KH}_2\text{PO}_4$ .

For TYA-agar:

Add 15 g of agar per liter.

For TYA-top agar:

Add 7.5 g of agar per liter.

Autoclave.

Last modifications: 2011-02-25